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Undecalcified bone alizarin Red stain report**I. Experimental equipment and reagent**

Equipment

Brand name	Manufactures	Model
Oven	Tianjin Labotery instrument equipment Co. Ltd	GFL-230
Microslide	Wanwu	
Optical Microscopy	NIKON	Nikon Eclipse E100
Imaging system	NIKON	Nikon DS-U3

Reagent

Reagent	Manufactures	Cat
Ethanol absolute	Sinopharm Chemical Reagent Co.,Ltd	100092183
Xylenes	Sinopharm Chemical Reagent Co.,Ltd	100023418
Alizarin Red staining solution	Wanwu	G1038
Neutral balsam	Sinopharm Chemical Reagent Co.,Ltd	10004160
Ethylene glycol ethyl ether acetate	Macklin	E808814-2.5L

II. Experiment procedure**1. Remove the embedding agent and rehydration**

Put sections into ethylene glycol ethyl ether acetate I 6 h at 37 °C and ethylene ether acetate II overnight at 37 °C. Then put them into ethylene glycol ethyl ether acetate III 10-15 minutes at room temperature and ethylene glycol ethyl ether acetate IV 10-15 minutes at room temperature. The sections were rehydrated in 100%I% - 100%II - 95% - 90% - 80% alcohol. Each step takes anywhere for 10 minutes. Finally, rinse with running water.

2. Alizarin Red staining

Put sections into alizarin red staining solution 5-10 minutes and wash with running water. Then place in the oven to drying.

3. Transparent and mount

Put sections into three cylinders of anhydrous ethanol for 5 min and two cylinders of xylene for 5 min. Finally, mount the sections with neutral balsam.

4. Microscope observation and collection and analysis images.**III. Result determination**

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Hefei WANWU technology CO.,LTD

Calcium salt deposits are crimson with a reddish background or near colorless.