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Climbing film, dropping film, experimental report

1. Experimental equipment and reagents

1.1 Experimental equipment

Name	Manufactor	Model
Six-well plate		
Dissecting forceps		
Pasteur pipette	Wanwu	WGXG0005
Bottle washing		
oven	Tianjin Lai Bo Rui Instrument Equipment Co., Ltd.	GFL-230
Glass slide	Wanwu	
Upright optical microscope	Nikon Japan	Nikon Eclipse E100
Imaging system	Nikon Japan	Nikon DS-U3

1.2 Main experimental reagents

Reagent name	Manufactor	Article number
Membrane Breaker	Wanwu	G1204
Fixative	Wanwu	G1101
Neutral gum	Sinopharm Group Chemical Reagent Co., Ltd.	10004160

2. Experimental steps

Cell slide staining:

2.1 Pre-treatment of climbing piece:

- ① Fixing of climbing piece: add 3ml of fixing solution to the hole and fix it for 15min. Absorb the fixing solution in the hole and wash it 3 times with pure water.
- ② Breaking film of climbing piece: add 3ml of broken piece to the hole Membrane liquid, membrane breakage for 15 min. Absorb membrane breakage liquid in the pores and wash with pure water 3 times.

2.2 Climbing slide staining: the subsequent staining steps are the same as the paraffin section, the staining time is appropriately extended, and the paste or pipette is used to clean or replace the staining solution and suck up the previous staining solution.

2.3 Climbing film: add appropriate amount of alcohol or water (according to the type of dyeing) and pick up the film. The film rests on the edge of the hole. After absorbing the alcohol, it is

placed in a 65° C oven and dried for 15 minutes. Take it out of the oven. Slide the well plate, write the number of each well plate on the slide, after dropping a drop of neutral gum on the slide, use tweezers to pick up the slide with the corresponding number, and seal the cell face down.

2.4 Observed under microscope and took images.